

Spor-Klenz Concentrate Sterilant is a fast-acting liquid sporicide specifically formulated for sterilization and disinfection of environmental, hard non-porous surfaces. Spor-Klenz Concentrate Sterilant contains the proven blend of peracetic acid, hydrogen peroxide and acetic acid to provide fast, effective microbial control of challenging spores such as *Bacillus subtilis* and *Clostridium sporogenes*. This formulation can be diluted with water, offering operational flexibility for use as a disinfectant or sanitizer. This product has an extensive documentation package to meet your validation objectives and is supported by the STERIS Technical Services team for individualized application and validation assistance.



## Spor-Klenz<sup>®</sup> Concentrate Cold Sterilant

Fast-acting sporicide in a versatile format

### FEATURES

- Concentrated, easy to dilute formulation
- Broad spectrum efficacy against bacteria, fungi, viruses and bacterial spores using AOAC protocols
- Excellent material compatibility
- Low toxicity profile
- 0.2 micron filtered and each lot USP sterility tested
- Manufactured in accordance with cGMP with complete lot traceability

### BENEFITS

- Allows use as a sporicide, disinfectant or sanitizer
- Reduces packaging waste and required storage space
- Confirms efficacy against many common microbes
- Allows use on most surfaces
- Supports worker safety
- Facilitates introduction into critical areas
- Supports quality assurance and quality control goals


### APPLICATIONS

Surface disinfection and microbial control of:

- Floors
- Walls
- Ceilings
- Countertops
- BSL hoods
- Carts and Other Equipment

Products may be applied by:

- Mopping
- Wiping
- Spraying

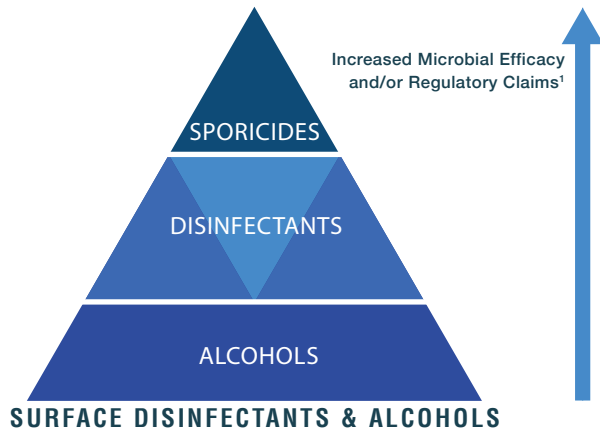

In maintaining microbial control of your facility, sporicides perform a critical function, destroying the toughest environmental contaminants. They should be used based on the frequency of spore results in your environmental data and should be used in rotation with one or more surface disinfectants. Spor-Klenz Concentrate Sterilant offers fast, broad spectrum efficacy with superior material compatibility. Our Technical Service group can help customize your contamination control program to take advantage of this excellent product.

<sup>1</sup>Products that fall into the categories at the bottom of the pyramid are most frequently used and are generally not sporicidal. Progression up the pyramid indicates stronger performance overall and a broader spectrum of claims.

## ORDERING INFORMATION

### Spor-Klenz Concentrate Cold Sterilant

Filtered for consistent quality

		United States
32 oz (946 mL) bottle	4 per case	652026